# Nano PCR <u>Purification Kit</u>



# **User Guide**



# ADVANCED MICRODEVICES PVT. LTD.

21, Industrial Area, Ambala Cantt - 133006 (INDIA)
Tel: 0171-2699290, 2699471, 2699274, Fax: +91-171-2699221, 2699008
Email: info@mdimembrane.com, Website: www.mdimembrane.com

# Table of Contents

		Page
1.	Introduction	3
2.	Applications	3
3.	Storage Conditions	3
4.	Quality Assurance	4
5.	Safety Information	4
6.	Lot Release Criteria	4
7.	Technical Support	4
8.	Kit contents	5
9.	Specifications	5
10.	Yields	5
11.	Procedure	6
12.	Principle	7
13.	Important Points to be Considered	8
14.	Protocol	9
15.	Trouble Shooting Guide	10
16.	Product Use Limitations	11
17.	Product Warranty and Satisfaction Guarantee	11
18.	Ordering Information	12

#### 1. Introduction

mdi Nano PCR Purification Kit is a fast, economical and easy way to obtain highly concentrated, purified DNA fragments (70bp-4kb) from PCR amplification reactions. The buffer system provided in the kit allows selective binding of DNA fragments from PCR reactions onto the spin column.

Washing is done with the help of provided wash buffer in order to remove primers, nucleotides, enzymes, mineral oil, salts and other impurities from DNA samples. Purified DNA is eluted in low salt buffer or water for a variety of downstream applications. The technology also does away with the cumbersome methodologies of phenol extraction (associated with slurries formation) as well as ethanol precipitation (associated with anion exchange based purification system) for desalting.

#### 2. Applications

- 1. Ligation
- 2. Transformation/Transduction/Transfection
- Automated Fluorescent Sequencing
- 4. Radioactive Sequencing
- 5. Restriction Digestion
- 6. Labeling
- 7. Cloning

## 3. Storage Conditions

As soon as you receive the kit, open it and store mdi Nano PCR spin columns at 2-8°C.

The remaining kit components can be stored at room temperature. The kit is stable for 1 year at given storage conditions. If precipitates appear in buffers, re-dissolve at 37°C for few minutes.

## 4. Quality Assurance

The mdi Nano PCR Purification kit is designed for various pre-determined specifications and user requirements such as yield, purity, ruggedness, shelf life and functional convenience.

These are produced through a well defined quality management system certified by Underwriters Laboratories, USA for ISO 9001: 2008 which ensures intra lot as well as lot to lot consistency.

#### 5. Safety Information

The buffers and the reagents may contain irritants, so wear lab coat, disposable gloves and protective goggles while working with the Nano PCR Purification Kit.

#### 6. Lot Release Criteria

Each lot of Nano PCR Purification Kit is tested against predetermined specifications to ensure consistent product quality.

### 7. Technical Support

At mdi, customers are our priority. We will share our experiences to assist you to overcome problems in general product usage as well as customize products for special applications. We will

- \* Stimulate problems, and suggest alternative methods to solve them.
- \* Make changes/improvements in our existing products/protocols.
- \* Develop special new products and systems especially to satisfy your needs.

We welcome your feedback to improve our products.

# 8. Kit Contents

Contents	Quantity			Storage Temperature
Spin Columns	50	250	1000	2-8°C
Buffer NPB	30ml	150ml	600ml	RT
Buffer PW	60ml	300ml	1200ml	RT
Buffer PE	15ml	75ml	300ml	RT
Collection Tubes	50	250	1000	RT
Hand Book	1	1	1	-
Certificate of Quality	1	1	1	-


# 9. Specifications

Capacity of column reservoir	800µl
Binding capacity of membrane (ds DNA)	5µg
Recovery of DNA	80%
Recovered DNA fragment	(70 bp - 4 kb)
Minimum elution volume	5μΙ
Total eluate volume	4µl

# 10. Yields

Feed	Yield	Percentage Recovery
422.7 ng	351.3 ng	83%
107.2 ng	83.64 ng	77%
61.67 ng	45.68 ng	73%

### 11. mdi Nano PCR Purification Procedure

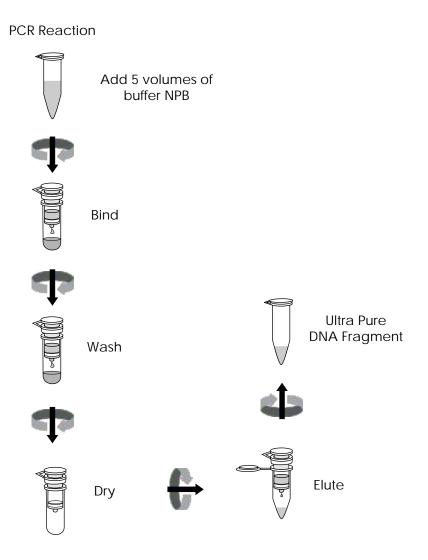

#### 12. Principle

The mdi Nano PCR purification kit allows the isolation of high ultra pure DNA fragments which involves:

- 1. Capturing PCR amplified DNA on spin column
- 2. Washing
- 3. Elution

#### 1. Capturing PCR amplified DNA on spin column

In order to facilitate adsorption of DNA fragments from PCR amplified reactions onto the spin column, optimum salt concentration and pH conditions are necessary which is achieved by addition of binding buffer 'NPB'.

#### 2. Washing

Subsequent to DNA binding onto the spin columns, unwanted components like primers, nucleotides, mineral oil, salts, and other impurities are washed away. Washing is done by buffer 'PW'.

#### 3. Elution

Salt concentration and pH of elution buffer is important for maximum elution efficiency, elution occurs at basic conditions and low salt concentration. Elution is done by buffer 'PE'.

#### 13 Important Points to be Considered

#### Optimization of Binding Buffers

The binding buffers posses appropriate salt concentration and pH to facilitate the efficient binding of single or double stranded DNA (70bp-4kb) from PCR reaction on spin column.

#### Centrifugation

\* All centrifugation steps should be carried out at room temperature at  $\geq$  10,000 rpm in a conventional table top microcentrifuge.

In case of choking of spin column, increase centrifugation time.

#### Washing

- \* To remove residual wash buffer spin the column for 1-2 minutes at ≥10,000 rpm.
- \* Flow through should be properly discarded before additional centrifugation step otherwise it will not be removed & may inhibit enzymatic reactions.

#### Elution

\* Elution buffer should be dispensed on to the center of column for maximum elution efficiency. If eluate volume is lower, then centrifuge for one minute extra at ≥ 10,000 rpm.

#### Yield

DNA Yield depends on following factors:

- 1. Volume of elution buffer
- 2. Point of application of buffer on the column
- 3. Incubation period of buffer on the column

#### 14. Protocol

Note: a) All steps should be carried out at room temperature.

- b) All centrifuge steps are at ≥10,000 rpm in a conventional table top micro centrifuge.
- 1. Add 5 volumes of buffer 'NPB' to 1 volume of PCR reaction and mix well. Users may or may not remove mineral oil or kerosene.
  - For example, add 250  $\mu$ l of buffer 'NPB' to 50  $\mu$ l PCR sample (not including oil).
- Place the spin column in 2ml collection tube in a suitable rack. Load the above sample solution in the center of the spin column, then centrifuge for 1min at ≥10,000 rpm and discard the flow through.
  - Note: Maximum volume of the column reservoir is 800 $\mu$ l. For sample volumes >800 $\mu$ l, simply load, balance and spin at  $\geq$ 10,000 rpm for 1min. Discard the flow through.
- 3. Place the spin column in the same collection tube. Wash the spin column with 750 µl of buffer 'PW' by giving a spin at ≥10,000 rpm for 1 minute. Discard the flow through.
- 4. Place the spin column in the same collection tube and centrifuge for 1-2 minutes at ≥10,000 rpm.
  - Important: This step is necessary to remove residual wash buffer. Residual wash buffer will not be completely removed unless the flow through is discarded before this additional centrifugation. Residual wash buffer may inhibit subsequent enzymatic reactions.
- 5. Place the spin column in a fresh 1.5ml microcentrifuge tube (not provided), elute DNA by dispensing 5µl of elution buffer 'PE' directly onto the center of the column membrane to maximize yields, allow it to stand for 1 minute and centrifuge for 1 minute at ≥10,000 rpm.

Note: The average eluate volume is  $4\mu l$  from  $5\mu l$  elution buffer. Centrifuge for one minute extra at  $\geq 10,000$  rpm if the eluate volume is lower.

#### 15. Trouble Shooting Guide

This section highlights some of the commonly faced problems, the probable reasons and suggestions to solve them.

Problem #1: Low or no recovery

Reason Elution buffer not properly

dispensed

Solution The elution buffer must be added

to the center of the membrane to ensure that the entire membrane area is covered by it and no elution buffer lies wasted sticking to the sides of the spin column and incubate 2-3 mins more

before centrifugation.

Problem #2: Poor quality of recovered DNA.

Reason Residual wash buffer in the spin

column.

Solution Incomplete drying results in wash

buffer in the eluate which is difficult to load in gel and also results in reduced yield and performance in down stream

applications.

Dry for one minute extra by centrifuging at ≥10,000 rpm to completely remove the wash

buffer.

#### 16. Product Use Limitations

mdi kits are developed and manufactured for research purpose only. The products are not recommended to be used for human, diagnostics or drug purposes for which these should be cleared by the concerned regulatory bodies in the country of use.

#### 17. Product Warranty and Satisfaction Guarantee

All mdi products are guaranteed and are backed by our

- a. Technical expertise and experience of over 30 years.
- b. Special mdi process for consistency and repeatability.
- c. Strict quality control and quality assurance regimen.
- d. Certificate of analysis accompanied with each product.

mdi provides an unconditional guarantee to replace the kit if it does not perform for any reasons other than misuse. However, the user needs to validate the performance of the kit for its specific use.

# 18. Ordering Information

To order please specify as below:

Туре		
Type	Code	
NPCK	NPCK	

XX

XX

Pack Size		
Pack Size	Code	
50	0050	
250	0250	
1000	1000	

0250

Example:

NPCK	XX	XX
111 011	,,,,	, , , ,



